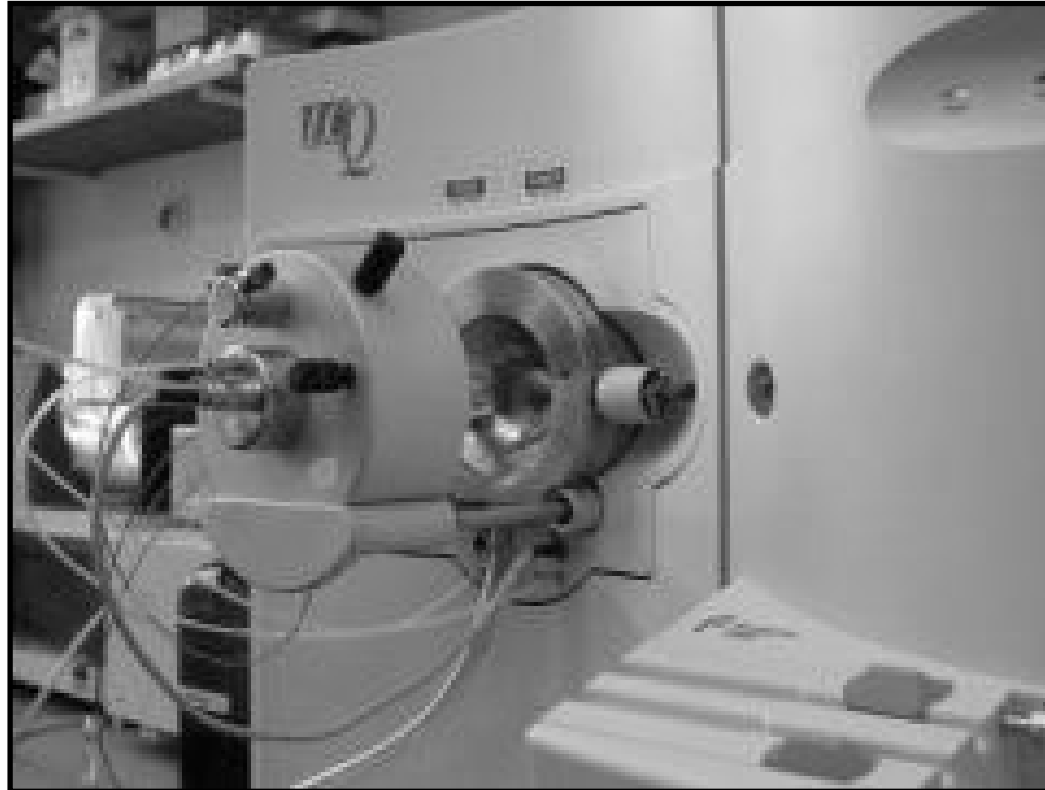
# Software Xcalibur 1.3

# LC-MS systém LCQ Classic

Thermo Finnigan

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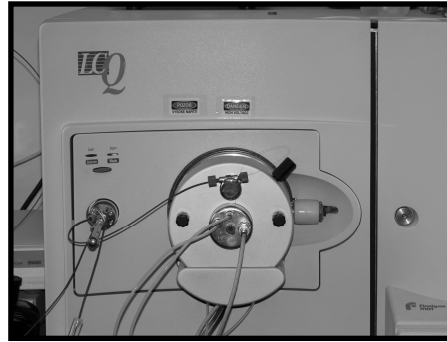
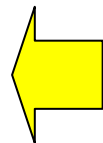
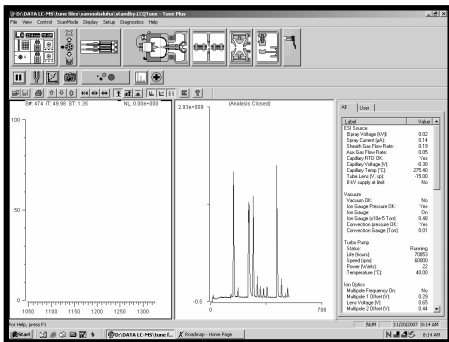
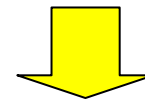
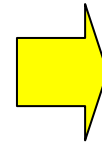
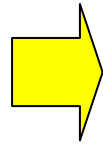
The screenshot shows the Tune Plus software interface in standby mode. The interface includes a menu bar (File, View, Control, ScanMode, Display, Setup, Diagnostics, Help), a toolbar with various icons, and a main display area with two plots. The left plot shows a signal trace with parameters S#: 474, IT: 49.98, ST: 1.35, NL: 0.00e+000. The right plot shows a mass spectrum with a peak at 2.93e+009. A status bar at the bottom shows 'Running' and '70853'.

Red arrows point from the following text boxes to specific icons in the toolbar:

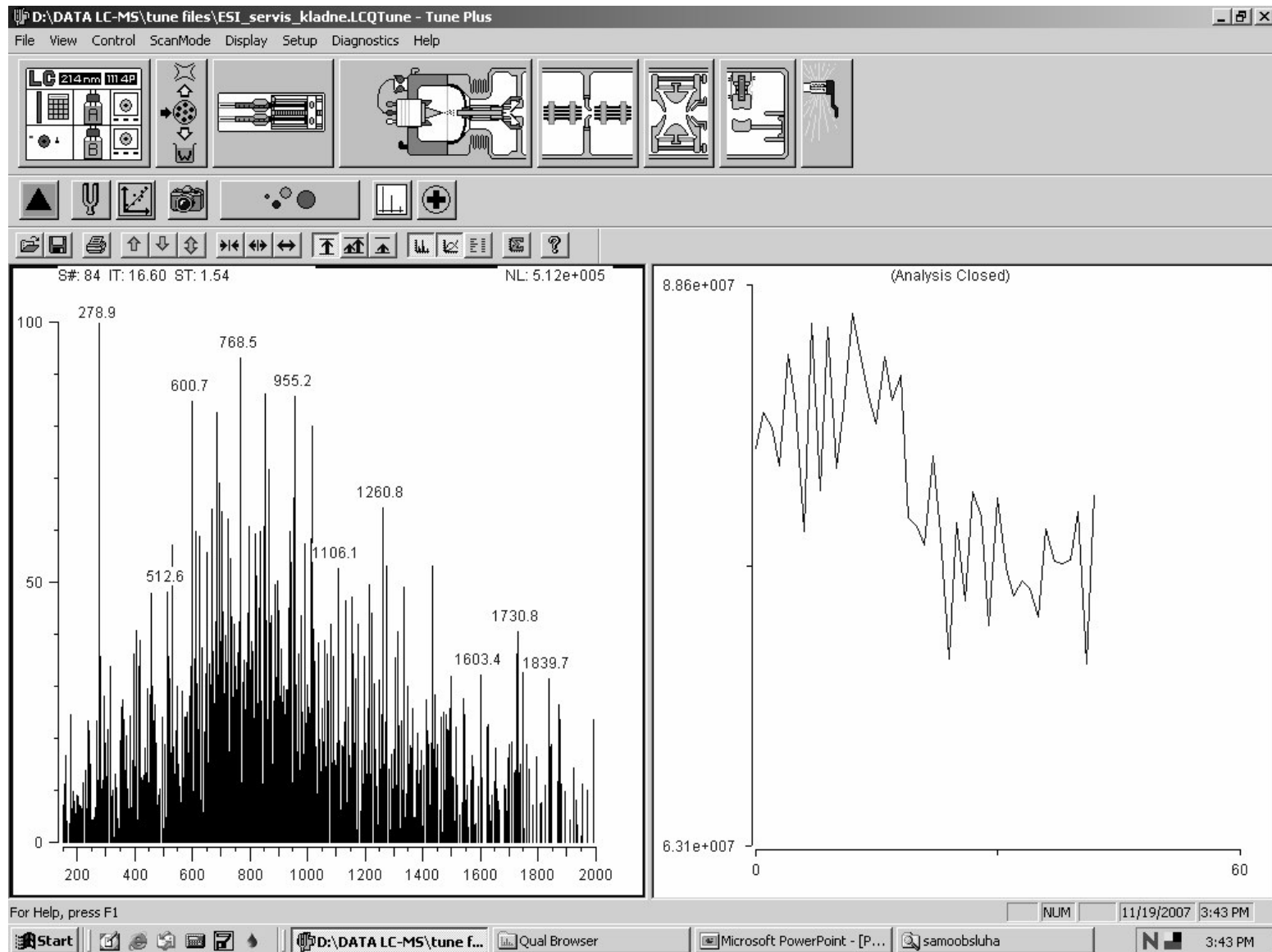
- Spouštění skenování (Start scanning) - points to the pause icon.
- Spouštění sběru dat (Start data acquisition) - points to the spray nozzle icon.
- Spouštění sledování signálu (Start signal monitoring) - points to the graph icon.
- Nastavení rozsahu a typu skenu (Set scan range and type) - points to the zoom icon.
- Typ dat : centroid/ profile (Data type: centroid/profile) - points to the data type icon.
- Volba polarity (nastaveno v ladící metodě) (Polarity selection (set in tuning method)) - points to the plus sign icon.

Label	Value
ESI Source	
Spray Voltage (kV):	0.02
Spray Current (µA):	0.14
Sheath Gas Flow Rate:	0.19
Aux Gas Flow Rate:	0.05
Capillary RTD OK:	Yes
Capillary Voltage (V):	-0.30
Capillary Temp (°C):	275.40
Tube Lens (V, sp):	-15.00
8 kV supply at limit:	No
Vacuum	
Vacuum OK:	No
Ion Gauge Pressure OK:	Yes
Ion Gauge:	On
Ion Gauge (x10e-5 Torr):	0.48
Convectron pressure OK:	Yes
Convectron Gauge (Torr):	0.01
Running	70853
	60000
	22
	40.00
Ion Optics	
Multipole Frequency On:	No
Multipole 1 Offset (V):	0.29
Lens Voltage (V):	0.65
Multipole 2 Offset (V):	0.44

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The screenshot displays the Tune Plus software interface. The menu bar includes File, View, Control, ScanMode, Display, Setup, Diagnostics, and Help. The File menu is open, showing options like Open..., Save, and Save As... with keyboard shortcuts. A list of tune files is visible, including '1 ESI\_servis\_kladne' and '5 D:\DATA LC-MS\...\draca\_APCI'. The main window shows a mass spectrum plot with peaks at 1866.8 and 657.8. A callout box highlights the 'Open' option in the File menu and the path 'D:\DATA LC-MS\tune files\samoobsluha'. The right panel displays instrument parameters such as ESI Source, Vacuum, Turbo Pump, and Ion Optics.

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Label	Value
ESI Source	
ISpray Voltage (kV):	4.53
Spray Current (µA):	0.39
Sheath Gas Flow Rate:	57.96
Aux Gas Flow Rate:	19.60
Capillary RTD OK:	Yes
Capillary Voltage (V):	11.42
Capillary Temp (°C):	275.40
Tube Lens (V, sp):	-15.00
8 kV supply at limit:	No
Vacuum	
Vacuum OK:	Yes
Ion Gauge Pressure OK:	Yes
Ion Gauge:	On
Ion Gauge (x10e-5 Torr):	0.90
Convector pressure OK:	Yes
Convector Gauge (Torr):	0.84
Turbo Pump	
Status:	Running
Life (hours):	70836
Speed (rpm):	60000
Power (Watts):	56
Temperature (°C):	45.00
Ion Optics	
Multipole Frequency On:	Yes
Multipole 1 Offset (V):	-2.71
Lens Voltage (V):	-15.15
Multipole 2 Offset (V):	-6.37

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The screenshot displays the Tune Plus software interface. The main window shows a chromatogram with several peaks labeled with their retention times: 278.8, 491.1, 651.3, 721.3, 1046.4, 1234.9, 1527.6, 1657.2, and 1879.2. The x-axis represents time in minutes, ranging from 0 to 2000. The y-axis represents intensity, ranging from 0 to 100. A red arrow points from the 'Inlet Direct Control' dialog box to the chromatogram. The dialog box is titled 'Inlet Direct Control' and contains the following settings:

Parameter	Value	Unit
A (%)	90	n/a
B (%)	10	n/a
C (%)	0	n/a
D (%)	0	n/a
Flow (ml/min)	0.20	n/a

The dialog box also includes buttons for 'Start', 'Stop', 'Help', 'Close', and 'Help'. The software interface includes a menu bar (File, View, Control, ScanMode, Display, Setup, Diagnostics, Help) and a toolbar with various icons for instrument control and data analysis. The status bar at the bottom shows the current date and time: 11/19/2007 3:35 PM.

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The screenshot displays the LCQ Tune Plus software interface. The main window shows a mass spectrum plot with the following parameters: S#: 1861, IT: 18.78, ST: 1.54. The x-axis represents m/z from 200 to 2000, and the y-axis represents relative intensity from 0 to 100. Key peaks are labeled with their m/z values: 358.3, 278.8, 702.1, 751.2, 958.6, and 1094.3. A 'Tune' dialog box is open in the foreground, with the 'Automatic' tab selected. The 'What to Graph' section has three radio buttons: 'TIC' (selected), 'ESI Spray Current', and 'Mass(es) (m/z)'. Below this, there are four checkboxes for mass values: 1 [ ] 195.10, 2 [ ] 524.30, 3 [ ] 1222.00, and 4 [ ] 1822.00. The 'Status' section is empty. At the bottom of the dialog are 'Start', 'Cancel', 'Print', and 'Help' buttons. A red arrow points from the 'Tune' dialog box to the 'ESI Spray Current' radio button in the software's toolbar. The Windows taskbar at the bottom shows the Start button, a file explorer window for 'D:\DATA LC-MS\tune f...', and a Microsoft PowerPoint window. The system tray shows the date and time as 11/19/2007 3:36 PM.

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The screenshot displays the Tune Plus software interface. The main window shows a chromatogram with several peaks labeled with their retention times: 278.7, 415.0, 544.2, 7, 1090.5, 1312.9, 1525.3, 1740.9, and 1997.3. The y-axis represents intensity, ranging from 0 to 100. The x-axis represents time in minutes, ranging from 200 to 2000. A red arrow points from the 'Acquire Data' dialog box to the 'Start' button in the main window's toolbar.

The 'Acquire Data' dialog box is open, showing the following fields and options:

- Filename: D:\DATA LC-MS\data\samoobsluha\inicialy\_vzorek.RAW
- Sample Name: popis vzorku
- Comment: +ESI, MeOH; 275C; sample in MeOH
- Acquire Time:  Continuously,  Scans (10),  Minutes (1.00)
- Acquisition Status: State: Idle, Time (min): 0.00
- Buttons: Start, Pause, View..., OK, Cancel, Help

The status bar at the bottom of the window shows the date and time: 11/19/2007 3:40 PM. The taskbar at the bottom of the screen shows the Start button and several open applications, including 'D:\DATA LC-MS\tune f...', 'Qual Browser', 'Microsoft PowerPoint - [P...', and 'samoobsluha'.

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D:\DATA LC-MS\tune files\samoobsluha\ESI\_servis\_kladne.LCQTune - Tune Plus

File View Control ScanMode Display Setup Diagnostics Help

LC 214nm M4P

**Define Scan**

S#: 318

Scan Description

Scan Mode:  MS  MS/MS  MSn

Scan Type:  Full  SIM  ZoomScan

MSn Power: [Slider]

TurboScan  Wideband Activation

MSn Settings

Parent Mass (m/z) Isolation Width (m/z) Normalized Collision Energy (%)

Scan Ranges (m/z)

From Mass To Mass

1 150.00 2000.00

Source Fragmentation

Turn On Collision Energy (V): 25.0

Apply OK Cancel Help

Nastavení módu a typu skenu

Nastavení rozsahu skenu

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NUM 11/19/2007 3:50 PM

Start D:\DATA LC-MS\tune f... Qual Browser Microsoft PowerPoint - [P... 3:50 PM


# MS/MS sken

The screenshot displays the 'Define Scan' dialog box in the Tune Plus software. The dialog is titled 'Define Scan' and contains several configuration sections:

- Scan Description:** Scan Mode:  MS/MS, Scan Type:  Full, MSn Power: 2.
- MSn Settings:** Parent Mass (m/z): 123.00, Isolation Width (m/z): 1.0, Normalized Collision Energy (%): 30.0. This section is highlighted with a red box and the text 'Nastavení parametrů MS/MS'.
- Scan Ranges (m/z):** From Mass: 50.00, To Mass: 200.00.
- Scan Time:** Number of Microscans: 3, Maximum Inject Time (ms): 200.00.
- Input Method:**  From/To,  Center/Width.
- Source Fragmentation:**  Turn On, Collision Energy (V): 25.0.

The background shows a mass spectrum plot with a peak at 278.9 m/z. The system status window on the right displays various parameters such as Voltage (kV), Gas Flow Rate, and Temperature.

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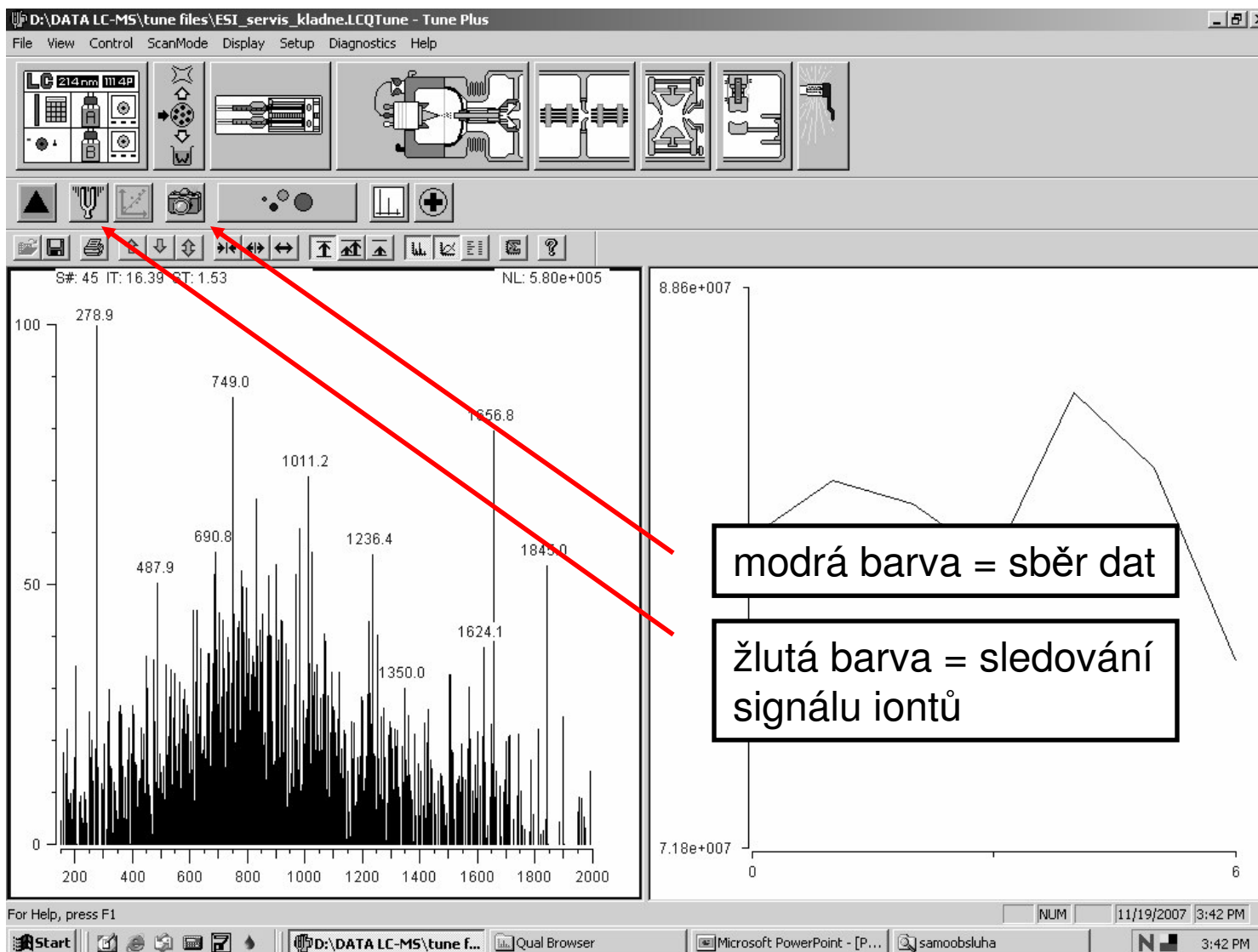# Okno programu Xcalibur

The screenshot displays the Xcalibur software interface. On the left, a sidebar titled "Acquisition Queue" lists the following components and their status:

- Run Manager
  - Ready To Download
  - Sequence:
  - Sample Name:
  - Working On:
  - Position:
  - Raw File:
  - Inst. Method:
- LCQ MS
  - Initializing
- TSP UV6000
  - Ready to Download
- TSP Autosampler
  - Not Connected
- TSP P4000
  - Ready to Download

The main workspace features a workflow diagram with the following steps:

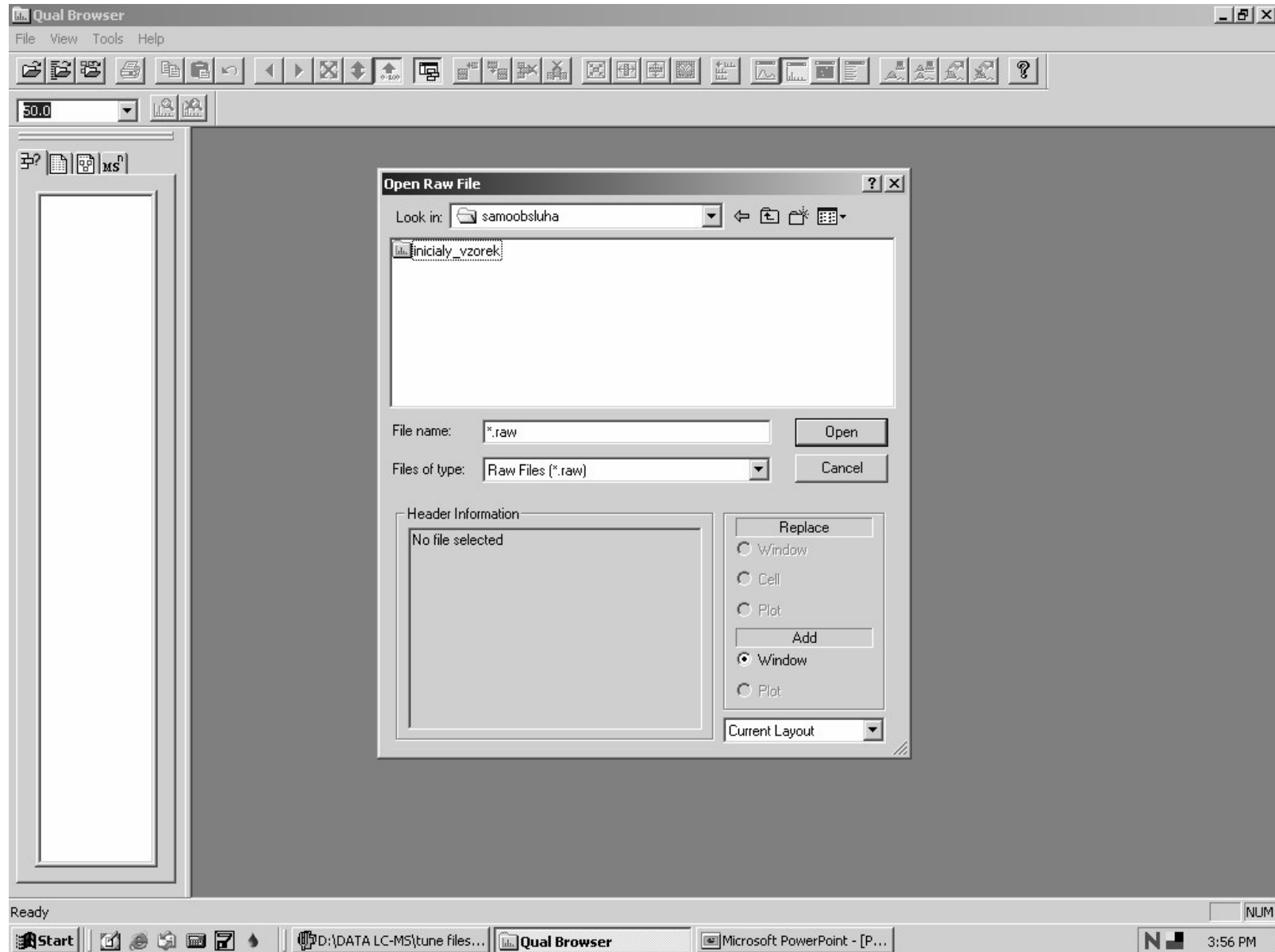
- Instrument Setup**: Represented by a keypad icon with "ON-LINE" text.
- Sequence Setup**: Represented by an icon of four vials.
- Processing Setup**: Represented by a gear icon.
- Results Review**: A large, semi-transparent watermark text.
- Qual Browser**: Represented by a molecular structure icon.
- Quan Browser**: Represented by a balance scale icon.
- Library Browser**: Represented by a chemical structure icon.

A red arrow points from a text box at the bottom left to the "Qual Browser" icon.

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
At the bottom of the window, the status bar shows "For Help, press F1" and the taskbar includes the Start button, a file explorer window for "D:\DATA LC-MS\tune files...", the "Roadmap - Home Page" window, and a "Microsoft PowerPoint - [m...]" window. The system tray shows the date "11/19/2007", time "4:32 PM", and the system tray icon.

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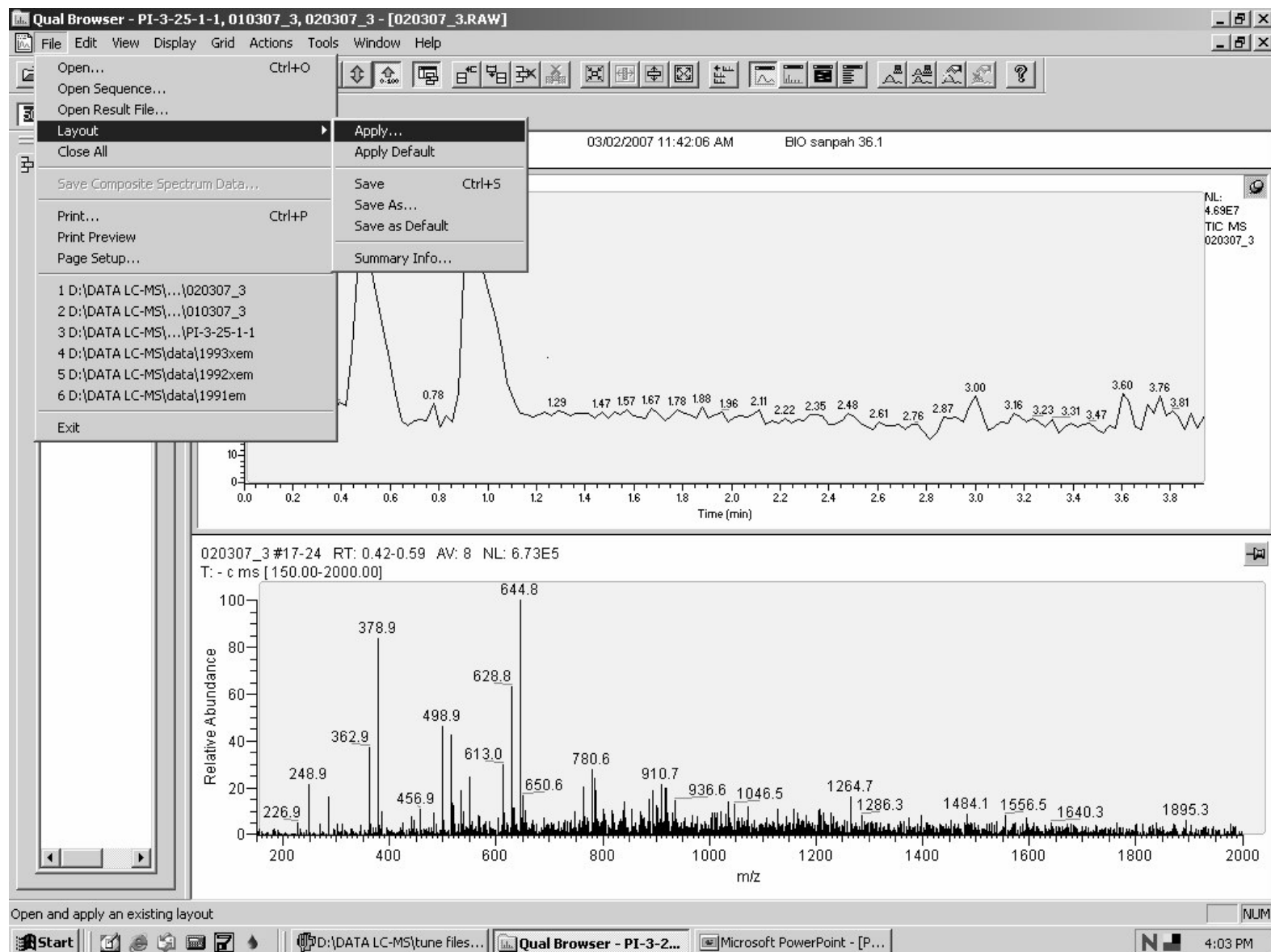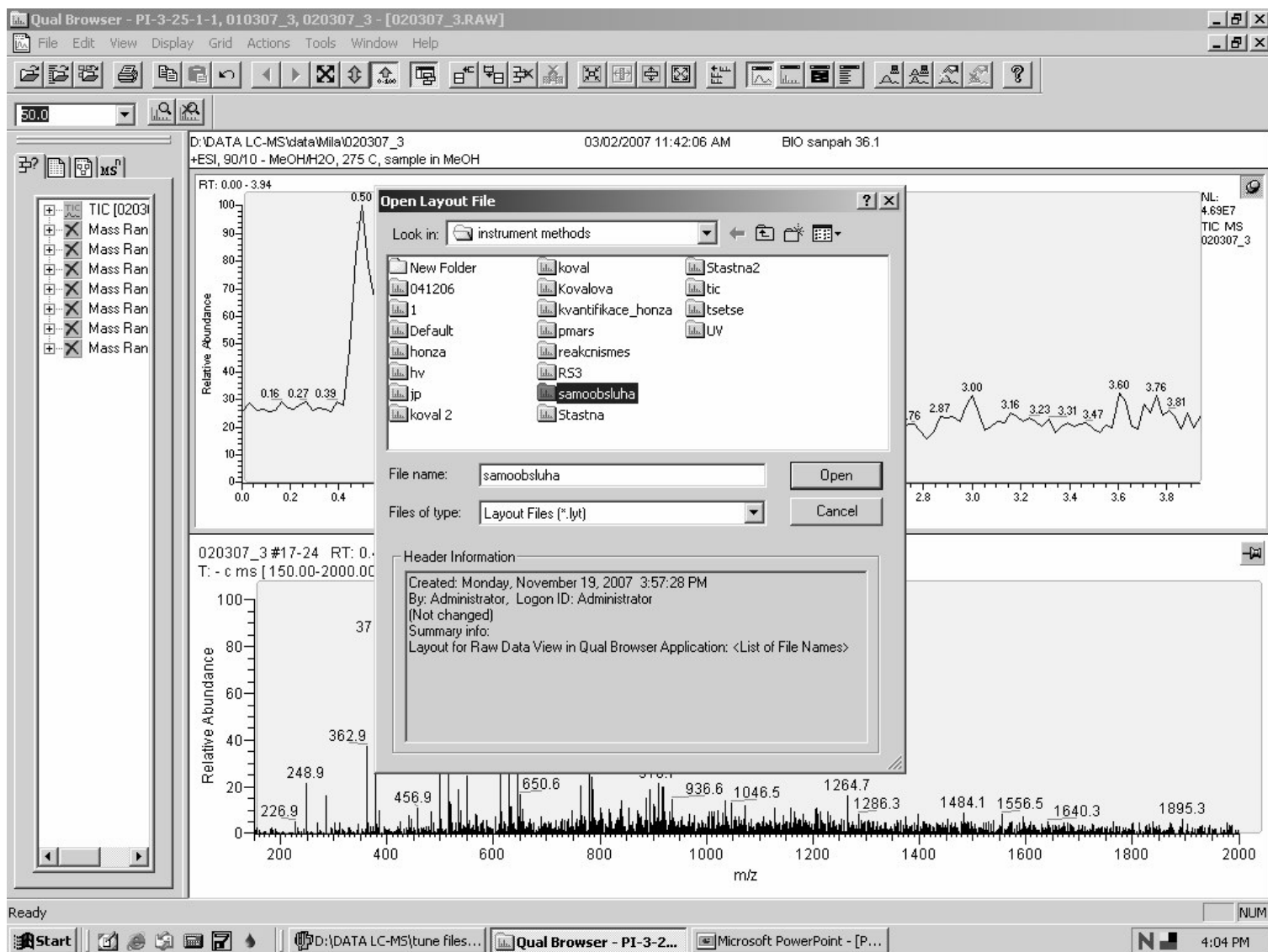
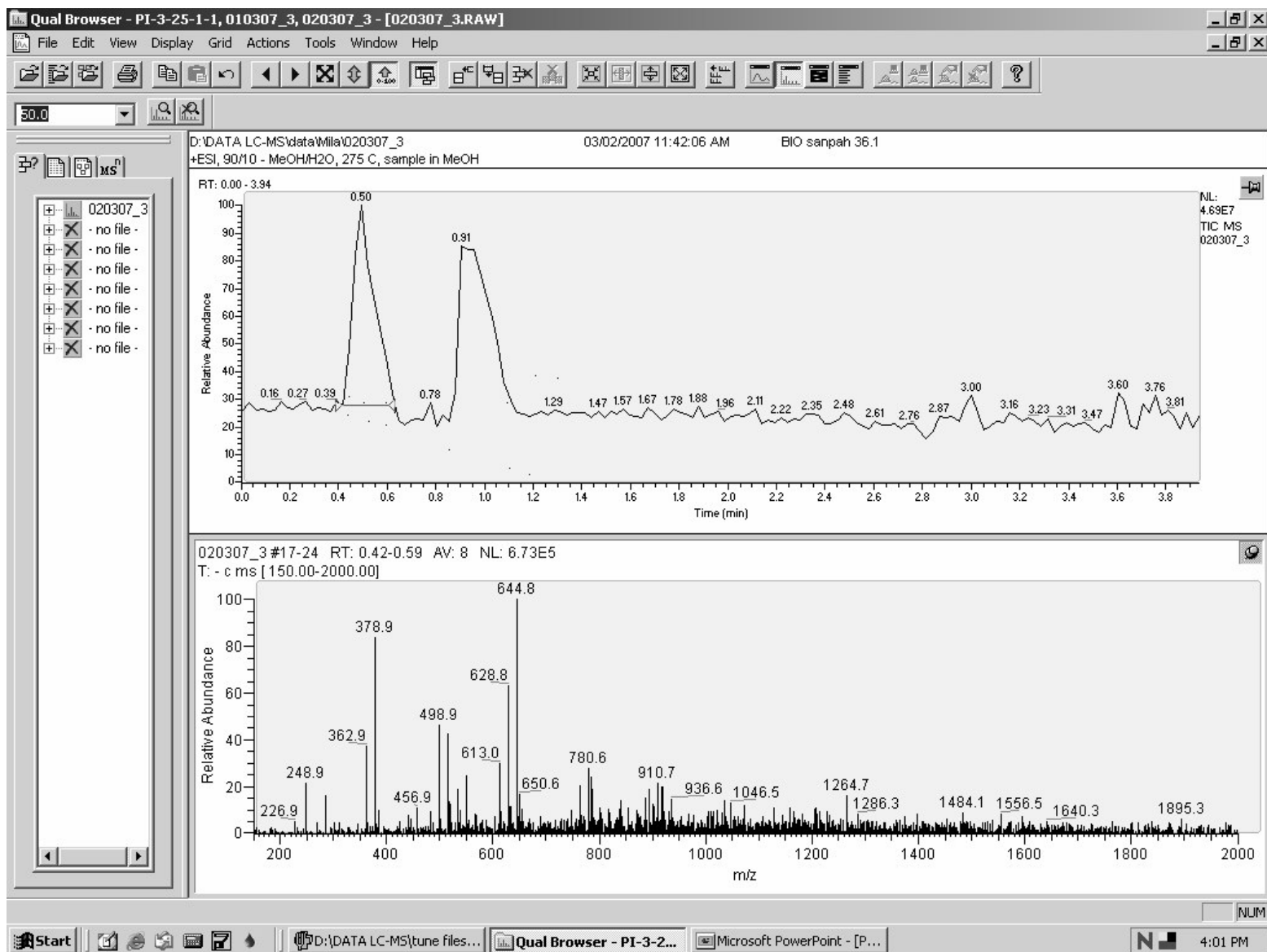
Qual Browser - PI-3-25-1-1, 010307\_3, 020307\_3 - [020307\_3.RAW]

File Edit View Display Grid Actions Tools Window Help

Subtract Spectra 1 Range  
Peak Detection 2 Ranges  
Peak Purity... Clear  
Library  
AutoFilter

D:\DATA  
+ESI, 90  
3  
sample in MeOH  
03/02/2007 11:42:06 AM BIO sanpach 36.1

RT: 0.00 - 3.94

Relative Abundance

Time (min)

NL: 4.69E7  
TIC MS  
020307\_3

020307\_3  
- no file -  
- no file -  
- no file -  
- no file -  
- no file -  
- no file -  
- no file -

020307\_3 #17-24 RT: 0.42-0.59 AV: 8 NL: 6.73E5  
T: - c ms [150.00-2000.00]

Relative Abundance

m/z

Subtract background scans, with 2 time ranges (usually before & after peak)

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