


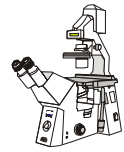
Manual

Cell Culture + Microscopy

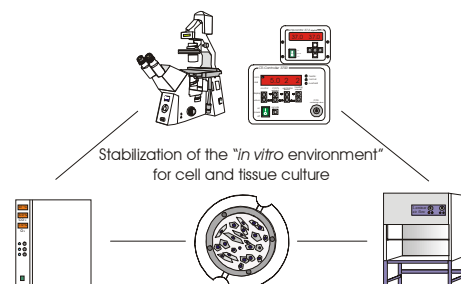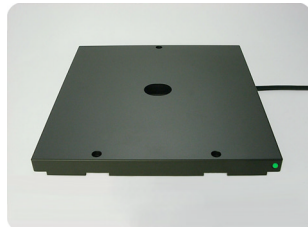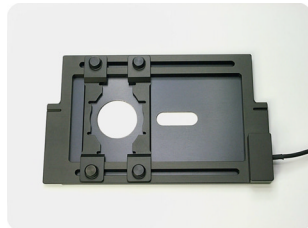Tempcontrol 37

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© 2001

Knowledge of this manual is required for the operation of the device. Would you therefore please make yourself familiar with the contents of this manual and pay special attention to hints concerning the safe operation of the device.

Design and specifications are subject to change without notice. The manual is not covered by an update service.

Unless expressly authorized, forwarding and duplication of this document, and the utilization and communication of its contents are not permitted. Violations will entail an obligation to pay compensation.

All rights reserved in the event of granting of patents or registration of a utility model.

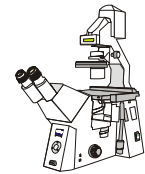
All product names mentioned herein may be the trademarks or registered trademarks of their respective companies and "TM" and "®" are not mentioned in each case in this manual.

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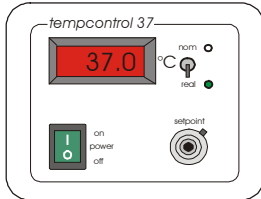
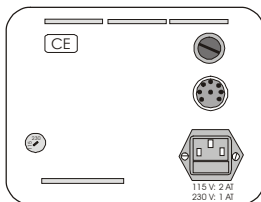
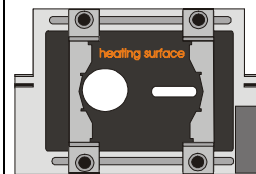
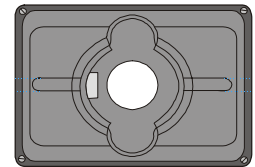
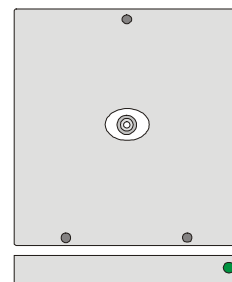
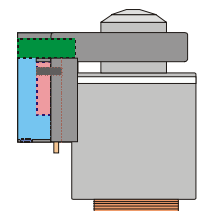
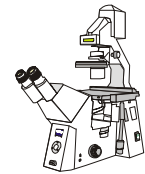
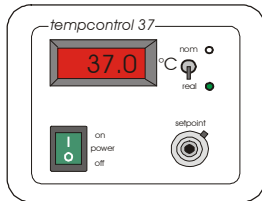
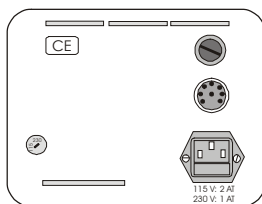
rear view

Description

- The temperature regulator Tempcontrol 37 (1-channel) is designed for heating and controlling of special microscope stages, plates etc. (see below).
- The device has one analogously controlled heating circuit (linear PID-regulator) and has a temperature control range from 3°C over room temperature to approx 60°C (adjusting range from 10°C up to 60°C).
- The real value is displayed on a digital multifunctional display (display in °C; resolution 0.1°C)
- This device is especially recommended for the control of Heatable Universal Mounting Frames.
- Controllable units (selection):

Heatable Universal Mounting Frames A-H, K-H und M-H

Heating Insert for stereo microscope Zeiss

Heatable Worktables

Heating Frame

CTI-Controller

Heating Microscope Stage

Heating and Cooling Frame

Heating Insert P and M 06-96

Objective Heater


Specifications

AC operating voltage	115 / 230 V, 50 ... 60 Hz
Rated power	90 VA max.
Fuse	115 V – 1.25 AT 230 V – 0.63 AT
Protective low voltage	24 V DC (output)
Control range	3°C over RT up to 60°C
Control tolerance	0.1°C
CE-Norm	class of protection 1
Dimensions (WxHxD)	175 x 138 x 200 (in mm)
Weight (kg)	3.8

Accessories (included)

Power cable # 0420.001

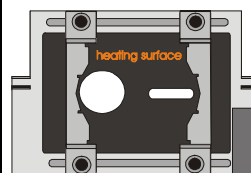Heatable Universal Mounting Frames M-H # 000000-1116-054


A-H # 000000-1116-055


K-H # 000000-1116-056

KH-L # 000000-1005-837

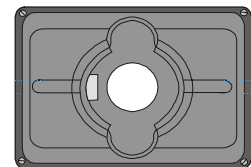Heating Insert P # 411861-9901-000

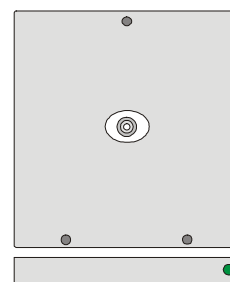Heatable Microscope Stage # 000000-1116-048


Objective Heater

000000-1116-073

000000-1116-075

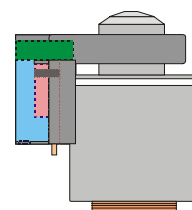000000-1116-076

Assembly:

- Ideally the device is placed on the right next to the microscope. The back side must have at least 10 cm free space to ensure sufficient cooling.
- Plug-in the 8-pin connector of the unit to be controlled into the socket on the back and tighten it by clockwise rotation.

Operation

- Switch-on the Tempcontrol 37 with the main switch at the front.
- Set the nominal value by shifting the switch to "**nom**", unlock the potentiometer and turn it until the desired temperature is displayed.
- To display the real value shift the switch back to "**real**". Lock potentiometer.

Maintenance

- The device is service-free.