
HXP 120 V

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Compact Light Source

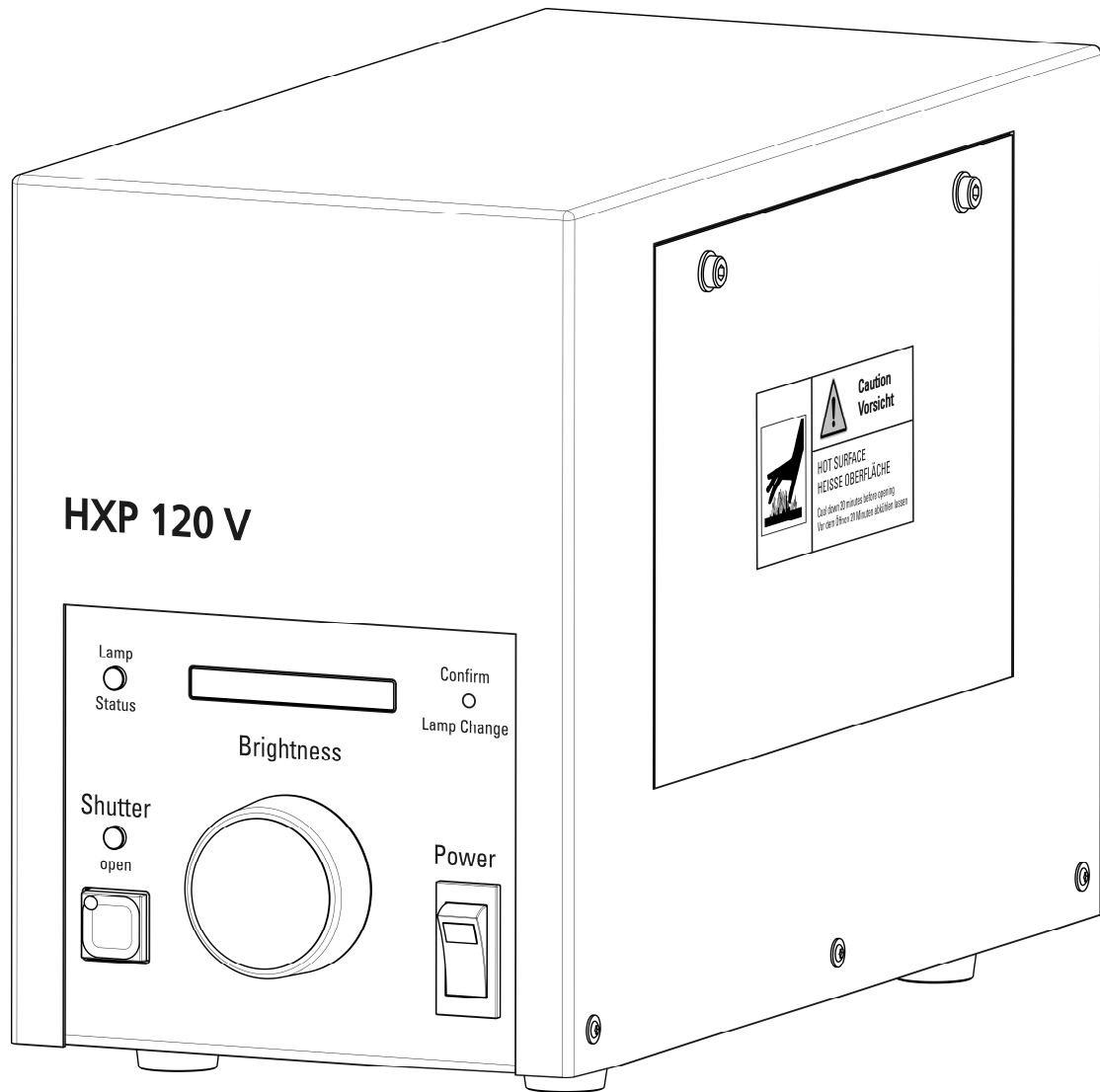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english

- © No part of this document may be transferred, copied by any method (print, photocopy, microfilm or other) or processed by electronic means unless with express permission. Any violation of this prohibition will be prosecuted.

The content of this document may change as design revisions of the product take place. These operating instructions are not updated..

Manufacturer/

editor:

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Stockholmer Straße 5

07747 Jena

Telephone no.: +49 (0) 3641 - 3530 0

Facsimile: +49 (0) 3641 - 35 30 70

Internet: www.lej.de

e-mail: info@lej.de

Closing date: 14. February 2013

Operating Instructions no.: 41 04 01-001-26A

Content

	Page
Copyright note	28
Content.....	29
Illustrations and tables.....	30
1	Fundamentals.....31
1.1	General information31
1.2	Product safety note32
1.3	Transport, storage and unpacking34
1.4	Terms of warranty34
2	Product description35
2.1	Intended purpose35
2.2	Product identification.....35
2.3	Scope of delivery35
2.4	Particular features of the product35
2.5	Connectible optical fibers36
3	Technical specifications37
4	Start-up41
4.1	Installation.....41
4.2	Installing and changing the lamp.....41
4.3	Protective filter44
4.4	Connecting the optical fiber44
4.5	Shutter remote and CAN-BUS45
4.6	Connecting the power cable46
5	Operation.....47
5.1	Controls47
5.2	Indicators48
5.3	Troubleshooting and remedies.....49
6	Service50
6.1	Maintenance and cleaning50
6.2	Spare parts and accessories50
6.3	Repairs51

Illustrations and tables

	Page
Figure 1	View of the Compact Light Source HXP 120 V 41
Figure 2	Device with housing cover removed; no lamp installed 42
Figure 3	Device with housing cover removed; lamp installed 43
Figure 4	Transmission characteristic of the protective filter 44
Figure 5	Rear of device with connections 45
Figure 6	View of controls of the Compact Light Source HXP 120 V 47
Figure 7	View of the indicators of the Compact Light Source HXP 120 V 48
Figure 8	Protective filter and filter glass 50
Table 1	Configuration "Trigger in" (Input) 39
Table 2	Configuration "Trigger out" (Output)..... 39
Table 3	Configuration "EXT" (Input)..... 39
Table 4	Faults and fault remedies 49
Table 5	Permitted fuses for use with the Compact Light Source HXP 120 V 51

Other

EC Declaration of conformity	52
------------------------------------	----

1. Fundamentals

1.1 General information

Read these operating instructions carefully before you connect and use the Compact Light Source HXP 120 V. The product must only be used by instructed persons. The user must also be familiar with the connection and operation of all components connected to the ballast. These operating instructions belong to the product and must be available for the life of the product. If the product is transferred to another user, the operating instructions must also be transferred.

The product was developed, manufactured and tested in accordance with norm EN 61010-1:2002 (IEC 1010-1) „Safety requirements for electrical equipment for measurement, control and laboratory use“ and in compliance with the applicable requirements of Annex 1 to Directive 2006/95/EG.

The Compact Light Source HXP 120 V complies with the requirements of EG directive 2004/108/EG. The norms and standards with which the product complies are listed in chapter 3 „Technical specifications“ in these operating instructions. Conformity with these norms and standards is documented by the CE label. The following hazard warning symbols have been used in these operating instructions:

**NOTE**

This symbol precedes information requiring your particular attention.

**NOTE**

Before the activity to which this note relates is carried out, disconnect all poles from the electric supply (pull mains plug).

**CAUTION**

This symbol warns of a hazard which can damage the product or any component (s) connected to it.

**CAUTION**

This symbol warns of a hazard which can cause injury to the user or other person.

**CAUTION**

Electric shock hazard!

**CAUTION**

Hazard due to high-energy light radiation. Do not look into the light beam. Use protective goggles.

**CAUTION**

Hazard due to hot parts. Allow to cool.

**CAUTION**

Explosion hazard Wear protective clothing.

1.2 Product safety note

The Compact Light Source HXP 120 V must only be used for the applications described in these operating instructions. Any other use is not covered by the manufacturer's warranty. If the device is used for a purpose for which it is not intended, all warranty and guarantee claims are voided. Further, the manufacturer's guarantee and warranty cannot be claimed if any of the components used with the Compact Light Source HXP 120 V is defective or unsuitable or if the product is serviced or repaired without authority.

If any safety defect or wrong function is found in the Compact Light Source HXP 120 V, the product must be switched off, all poles disconnected from the power supply and secured against use. Return the Compact Light Source HXP 120 V for repairs to the supplier or an authorized repair center.

The Compact Light Source HXP 120 V produces high-energy light some of which is not visible. There is risk of dazzling and blindness. Do not look in the optical fiber output of the device or the output of the optical fiber connected to the device.

Always make sure that the output of the connected optical fiber is firmly connected to the system which you want to illuminate before the compact light source is switched on.

The lamp used in the device becomes very warm in use (risk of burn) and has very high inside pressure when hot (explosion risk). Allow the device to cool for at least 20 minutes before you change the lamp.

The Compact Light Source HXP 120 V must only be connected to a functioning 2-pole and ground socket outlet. The product is only designed for use in a dry room. It is not permitted to use the product in a potentially explosive atmosphere.

Disconnect all poles of the device from the power line before connecting a component to the Compact Light Source HXP 120 V or starting service work which the user can do (lamp change; fuse change). If the fuses are changed, use only fuses of the type and rating specified in this document. Using makeshift fuse or rendering fuses ineffective (shorting the fuse holder, etc.) is forbidden.

The Compact Light Source HXP 120 V is designed for the voltage and frequency range of 100 ... 240 VAC, 50..60 Hz. In these voltage and frequency ranges, the product adjusts automatically to the power line voltage. Use of the compact light source outside the specified ranges can destroy the device and the connected components.

To avoid damage to the optical fiber connected to the device, the protective filter in the device (see „5 Operation“) must be free of defect (cracks, flaws, etc.). If the protective filter is not installed or is defective, the optical fiber which is connected to the device will be destroyed.

The lamp in the device contains highly toxic substances. At the end of the lifecycle, the lamp must be disposed of as prescribed. The lamp supplier provides the respective information (e.g., in the documentation supplied with the lamp). The documentation also contains supplier information about what to do if the lamp explodes.

1.3 Transport, storage and unpacking

To avoid damage to the lamp and the device, the lamp must be removed before the Compact Light Source HXP 120 V is transported.

The ambient conditions for transport and storage of the device specified in these operating instructions must be observed. It is safest to transport the Compact Light Source HXP 120 V in the original packaging. If damage is found when the Compact Light Source HXP 120 V is removed from the original packaging, inform the supplier or the forwarder without delay.

For unpacking, open the carton on top. The Compact Light Source HXP 120 V can be lifted from the carton together with the foam mold. Then the foam mold and the plastic bag can be removed from the device. Inspect the delivery of the compact light source (see „2,3 Scope of delivery“) for missing items. After unpacking, leave the Compact Light Source HXP 120 V for about 1 hour to adjust to room conditions before you start using it. This avoids damage from condensing water which forms when the temperature difference is high inside the device.

1.4 Terms of warranty

No repair or maintenance except as specified in these operating instructions must be carried out on the device. Repairs must only be carried out by the device supplier or an authorized repair center.

2 Product description

2.1 Intended purpose

The Compact Light Source HXP 120 V produces light of very high intensity and couples the light in the optical fiber, preferably liquid optical fiber with an active diameter of 3 mm. The lamp must be of the type specified in „3 Technical Specifications“. Any other use of the device is not intended.

2.2 Product identification

The nameplate of the device is located on the underside. Complete data for the identification of the device is printed on the nameplate.

2.3 Scope of delivery

The delivery of the Compact Light Source HXP 120 V contains the following components:

- Compact light source HXP 120 V
- Operating instructions
- Protective filter (integrated in the device)
- Liquid optical fiber (Lumatec Series 380, D=3, L=2000)
- BNC - cable
- Lamp OSRAM HXP-R120W/45C VIS
- Technical Information No. FO 5148

2.4 Particular features of the product

The Compact Light Source HXP 120 V has an automatic beam cover. The cover ensures that the light output of the device is closed if the optical fiber is pulled. This prevents that the user can be injured by the intensive light radiated by the device.

The Compact Light Source HXP 120 V has a built-in shutter. The shutter can be controlled by the „Shutter“ button at the device or remote controlled via several sockets.

The light source HXP 120 V has an integrated optical-mechanical dimmer unit. With this unit, the emitted brightness can be dimmed in 20 steps between 0 ... 100 % without changing other parameters of the light (e.g., light temperature). The dimmer unit can be set at a rotary switch at the device or remotely via CAN bus. An LED bar graph on the front of the device displays the brightness setting of the light source.

Built in the device is a filter holder with protective filter. The filter protects the optical fiber and the devices connected to it both from infrared radiation (heat protection function) and from ultraviolet radiation. The protection filter is exposed when the cover of the housing is removed and can be changed by the user if it is damaged.

2.5 Connectible optical fibers

Connected to the Compact Light Source HXP 120 V can preferably be liquid optical fibers with a light input of type „Storz, long“ and whose transmission characteristic matches the lamp used. The spectral light distribution of the lamp is specified in the data sheet of the lamp. The device is optimal for optical fibers with 3 mm active diameter.

Information on suitable optical fibers can be obtained from the supplier of the Compact Light Source HXP 120 V.

3 Technical specifications

Dimensions (width x height x depth)..... 145 mm x 210 mm x 315 mm
(incl. controls and connection elements)

Weight..... app. 5 kg

Ambient conditions:

Storage and transport (in original packaging):

Permitted ambient temperature.....-20 to 85°C

Permitted relative humidity 10 to 90 %, non-condensing

Operation:

Permitted ambient temperature.....0 to 40°C

Permitted relative humidity 10 to 90 %, non-condensing

Operating data:

Location Closed dry room

Altitude max. 2000 m ASL

Protection class I

Protection type IP20 (EN 60529)

Operating position standing upright

EMC:

Complies with..... EN 61326 -1 (2006-10)

The product complies with the requirements of EC directive 2004/108/EG

Electrical safety:

complies with.....EN 61010-1:2001

Pollution severity2

Overvoltage categoryII

Line power input:

Design Inlet connector for non-heating devices IEC320 C13 (EN 60320 C13)

Line power voltage 100 to 240 VAC (±10%)


Line power frequency 50 to 60 Hz

Fuses 2x T2.5 A; switching capacity H

Power consumption max. 210 VA

Lamp:

Type Mercury short arc reflector lamp
 Model HXP-R120W/45C VIS
 Supplier OSRAM
 Consumption 120 W
 Mean life (supplier information) 2000 h
 Spectral distribution of the output power see chart
 (supplier information)

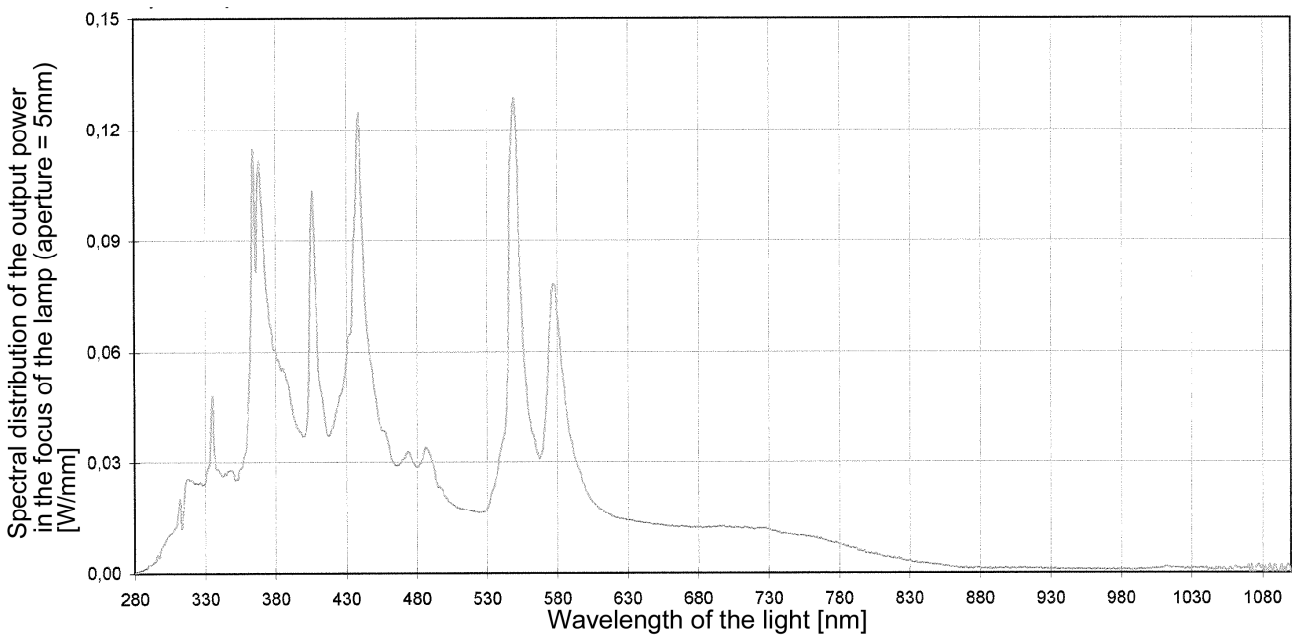Light output:

Suitable for adapter Type Storz, long
 Feature Self-closing if no optical fiber is plugged in
 optimal for active optical fiber diameter of 3 mm
 suitable, e.g., for Liquid optical fiber series 380 (product of LUMATEC)

Shutter:

Shutter frequency, max. 40 Hz
 Frequency limits > 45 Hz
 Maintenance-free operation (average) 2×10^7 cycles

Interfaces:

Trigger in:

Function Control of the integrated shutter
 level-triggered
 Design BNC connector 50 Ohm
 Level < + 0.4 V shutter close
 > + 3.0 V shutter open
 Max. input voltage 7 V
 Potential Screen on PE
 Configuration see table 1

Connection	Function
Inner conductor	Signal Trigger in
Screen	Signal Trigger in – GND

Table 1: Configuration "Trigger in" (Input)

Trigger out:

Function Shutter status output
 Design BNC connector 50 Ohm
 Level low-ohm \triangleq Shutter closed
 High-ohm \triangleq Shutter open
 Max. input voltage 20 V
 Max. current 3 mA
 Potential Screen on PE
 Configuration see table 2

Connection	Function
Inner conductor	Signal Trigger out
Screen	Signal Trigger out – GND

Table 2: Configuration "Trigger out" (Output)

EXT:

Function Control of the integrated shutter
 by button, flank-triggered
 Design BNC connector 50 Ohm
 Potential Screen on PE
 Configuration see table 3

Connection	Function
Inner conductor	Button Connection 1
Screen	Button Connection 2 – GND

Table 3: Configuration "EXT" (Input)

Protection functions:

Safety circuit

If the cover of the housing is opened, all components connected to the device (except power line input and power line switch) are disconnected from the power supply

Overtemperature protection

If overtemperature is detected, the lamp power supply is disconnected in the device. When the device has cooled, the power supply is established again automatically and the lamp fires.

Anti-dazzle

If no optical fiber is connected to the light output of the device, the light output is closed automatically and prevents the emission of light.

4 Start-up

4.1 Installation

 Set up the Compact Light Source HXP 120 V so that the front with the controls is freely accessible and visible. Install the compact light source on a safe and slip-proof surface. The ventilation slots at the sides and the rear of the device must not be covered, the free space in the area of the ventilation slots must at least be 150 mm. The cables and optical fibers connected to the Compact Light Source HXP120C must not create a risk.

 !

4.2 Installing and changing the lamp

 Caution: The lamp contains mercury. The handling and safety instructions supplied by the lamp manufacturer together with the lamp must be observed.

 Make sure that all poles are disconnected from the power line before the lamp is installed or a new lamp fitted in the Compact Light Source HXP 120 V. The Compact Light Source HXP 120 V must have cooled at least 20 minutes before the lamp can be installed. (Caution: The parts inside the device (lamp, etc.) take more time to cool than outside parts).


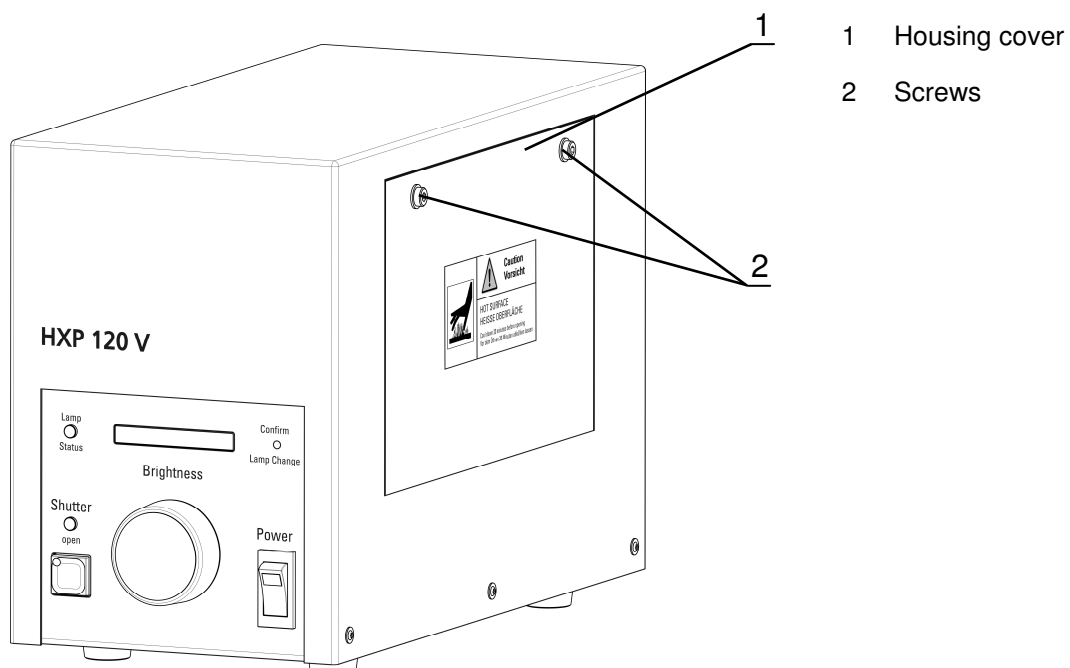
Figure 1: View of the Compact Light Source HXP 120 V

For installing the lamp, the housing cover (Figure 1[1]) of the compact light source must be removed. For this, remove the two screws (Figure 1[2]) with an Allen wrench WAF 3.0. If a lamp is fitted in the Compact Light Source HXP 120 V, that lamp must be removed before a new lamp can be installed. Remove the lamp in reverse order of the installation described here.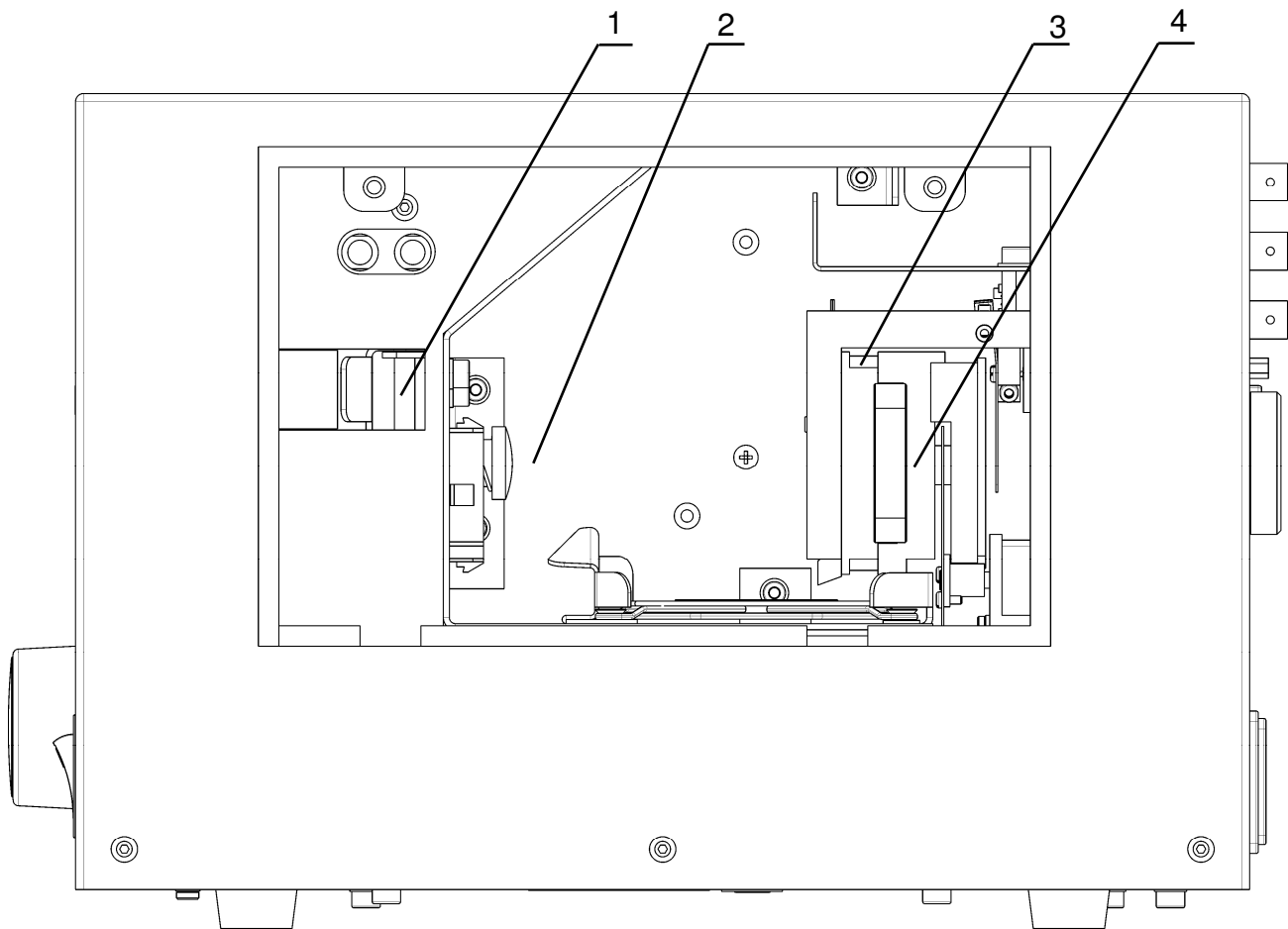


- 1 Lever
- 2 Pressure bolt
- 3 Lamp holder
- 4 Protective filter

Figure 2: Device with housing cover removed; no lamp installed

For installing the lamp, pull back the pressure bolt (Figure 2[2]) with the lever (Figure 2[1]) (move the lever towards the front panel) and install the lamp. Make sure that the groove in the contact face of the lamp engages the respective nose in the lamp holder (Figure 2[3]). If the lamp makes full contact with the lamp holder (Figure 2[3]), the lever (Figure 2[1]) can

slowly be moved towards the lamp. Make sure that the connecting cables of the lamp (Figure 3[3]) do not contact the reflector or block the venting grate (Figure 3[5]). To ensure this, install the front connecting cable of the lamp (at the light exit side) behind the air baffle (Figure 3[4]). When the lamp is installed and the plug connection made, the housing cover can be closed.

For removing the lamp it is recommended to lightly force the two parts of the plug (Figure 3 [1.2]) apart with a suitable screwdriver because the connection is very tight. During this, press down the locking lever at the plug (Figure 3[2]). Apply the screwdriver in such a way that when it slips only the lamp-side plug is damaged (Figure 3[2]) and not the socket (Figure 3[1]) as the latter is replaced together with the lamp.

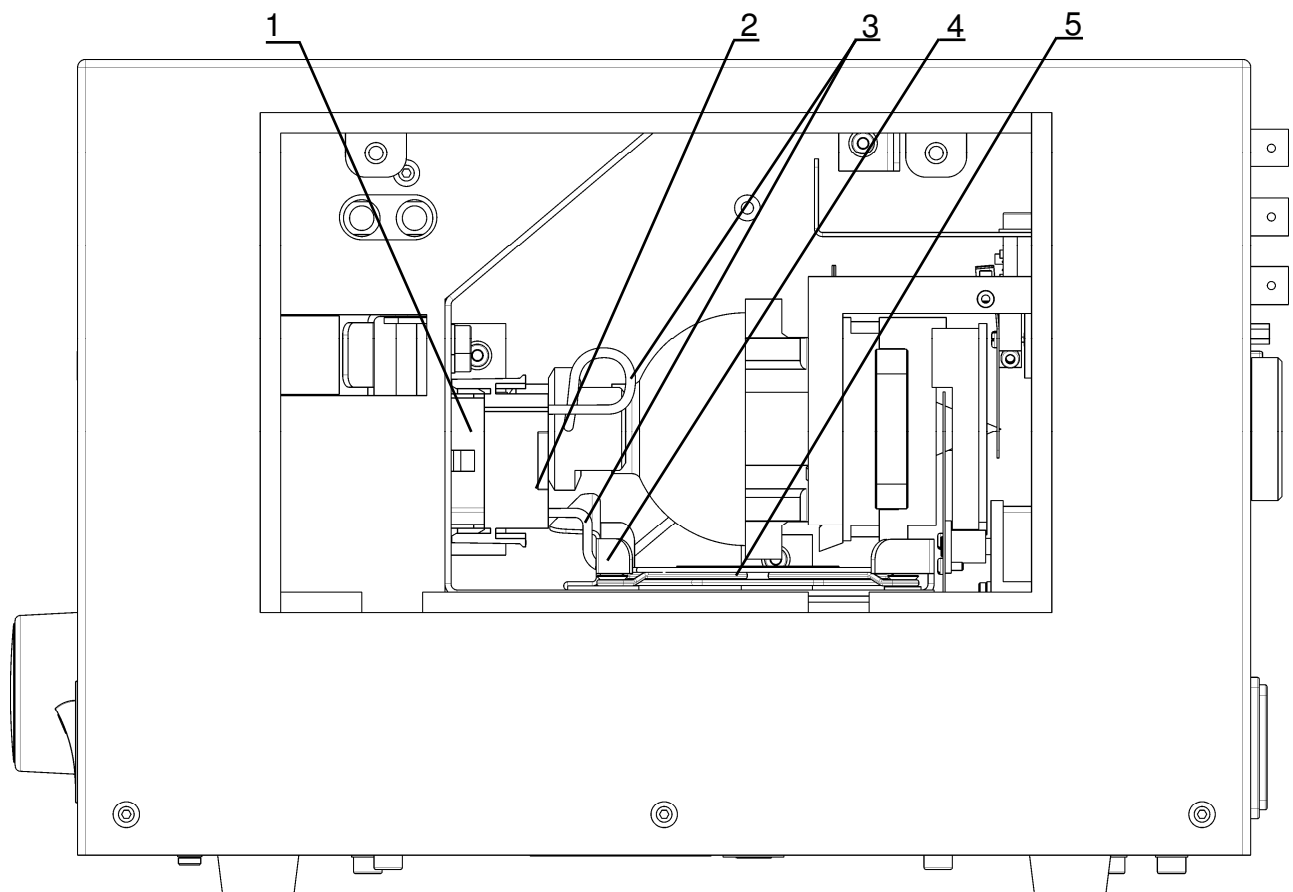- 1 Socket
- 2 Lamp-side plug

- 3 Connecting cables
- 4 Air baffle

- 5 Fan grate

Figure 3: Device with housing cover removed; lamp installed

4.3 Protective filter

To protect the optical fiber connected to the device and downstream devices from damage due to heat and ultraviolet radiation, a protective filter is integrated in the device (Figure 2 [4]). Figure 4 shows the transmission characteristic of the protective filter in the lamp. If the filter glass of the protective filter is damaged, it can be replaced by the user (see „6.2 Spare parts and accessories“). To replace the glass, the protective filter must be pulled out of the device. This can be done after removing the housing cover (Figure 1[1]) (see „4.2 Installing and changing the lamp“).

 Never use the device with a damaged or unsuitable protective filter or without a protective filter because the connected optical fiber and downstream devices can be damaged or destroyed.

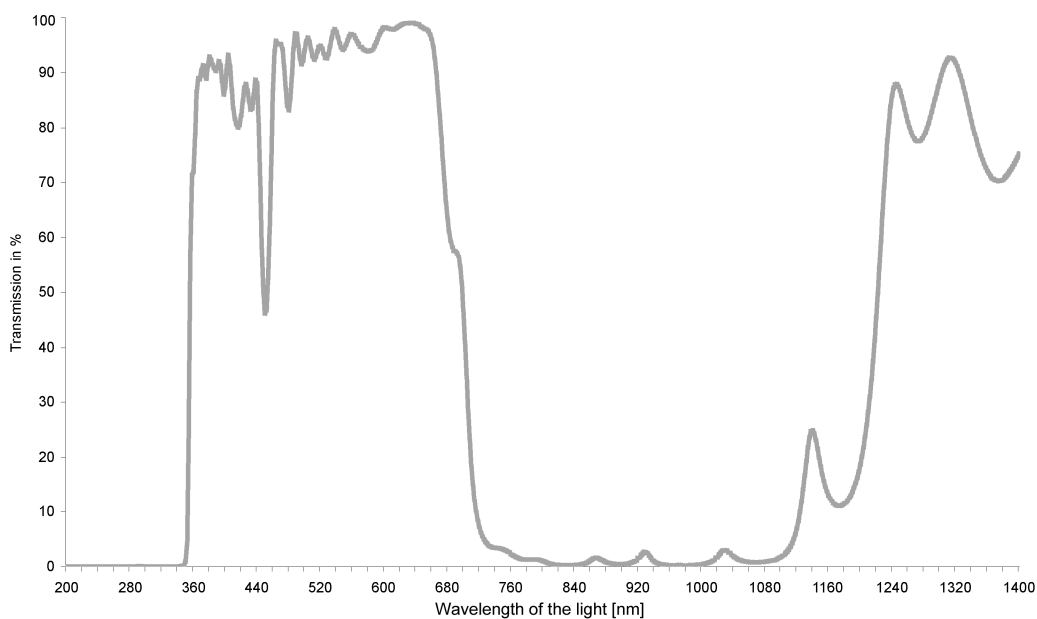
Figure 4: Transmission characteristic of the protective filter

4.4 Connecting the optical fiber

Always connect the output end of the optical fiber to the target system at first; this prevents hazard to which the user would be exposed due to the high-energy light emitted by the Compact Light Source HXP 120 V. Connect the input of the optical fiber to the light output at the rear of the compact light source (Figure 5[1]). The fiber should latch audibly. Only an optical fiber with light inlet of the type „Storz long“ can be used, otherwise the device can suffer damage and the user injury (by dazzling). Before you switch on the device, check again to make sure that the optical fiber is pushed in properly to the latching point.

4.5 Shutter control and CAN BUS

As alternative to the shutter on the device (Fig. 6[4]), the shutter in the device can be controlled via the connections „Trigger in“ (Fig. 5[2]), "EXT" (Fig. 5[4]) or the CAN BUS connection (Figure 5[5]).

It is possible to control the shutter via several connections at a time.

The „Trigger in“ input is only processed if the shutter was previously closed by another way (e.g., by pressing the shutter button). This input is level-triggered, which means that the shutter will not change position as long as the respective signal is active.

The „EXT“ input is provided for connection of a button. This input is flank-triggered, i.e., every time the button is pressed the shutter changes its position (like with the integrated shutter button).

Only devices which are equipped with the specific CAN BUS from Carl Zeiss (e.g., microscope) can control the shutter via the CAN BUS connection of the compact light source. This connection is designed as 15-pin HD SUB-D connector.

The connection should only be made or separated when the device is switched off.

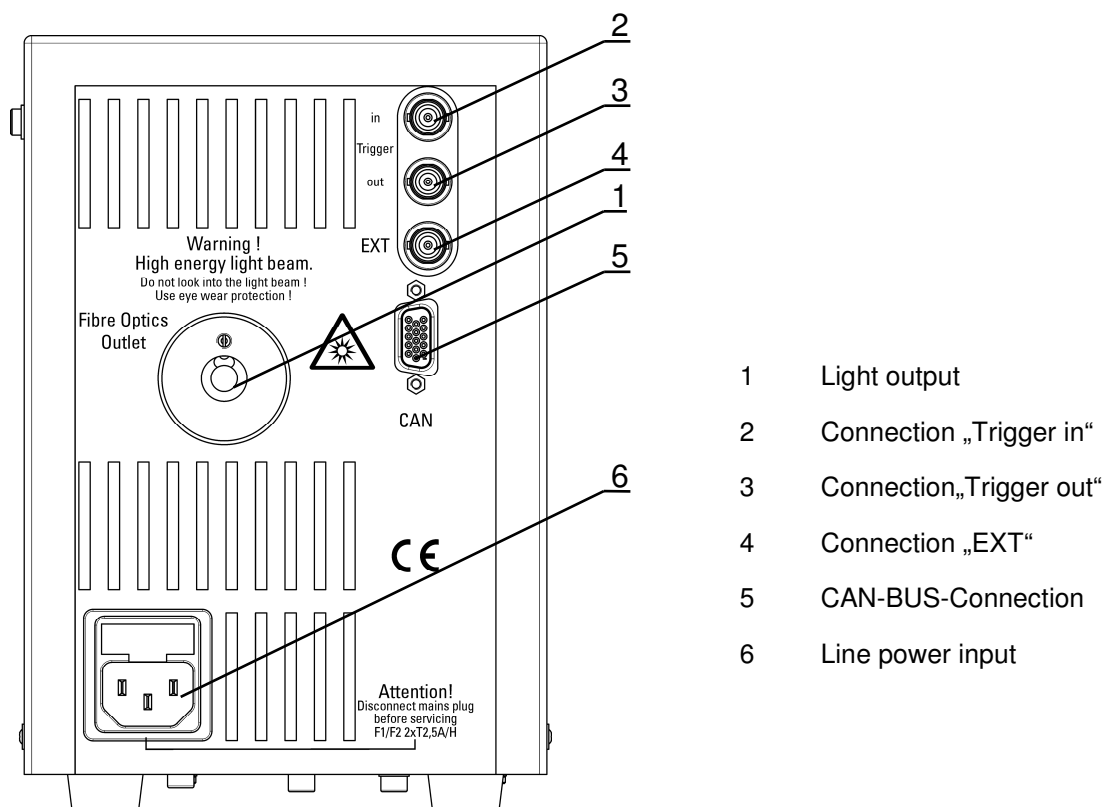
Figure 5: Rear of device with connections

4.6 Connecting the power cable

A power cable for non-heating appliances is used (conductor size: min. 3x0.75 mm²) with a socket outlet for non-heating appliances IEC320 C13 (EN 60320 C13) as output. The inlet plug of the connecting cable for non-heating appliances should conform to the site conditions. The protective conductor must be included in the connecting cable for non-heating appliances. Set the ballast mains switch to zero (Off) before connecting the mains cable.

At first, connect the mains cable to the Compact Light Source mains inlet (Figure 5 [6]) and then put the plug in a properly working ground socket outlet. The ground contact line must not be interrupted (e.g., by unsuitable extension cable).

5 Operation

5.1 Controls

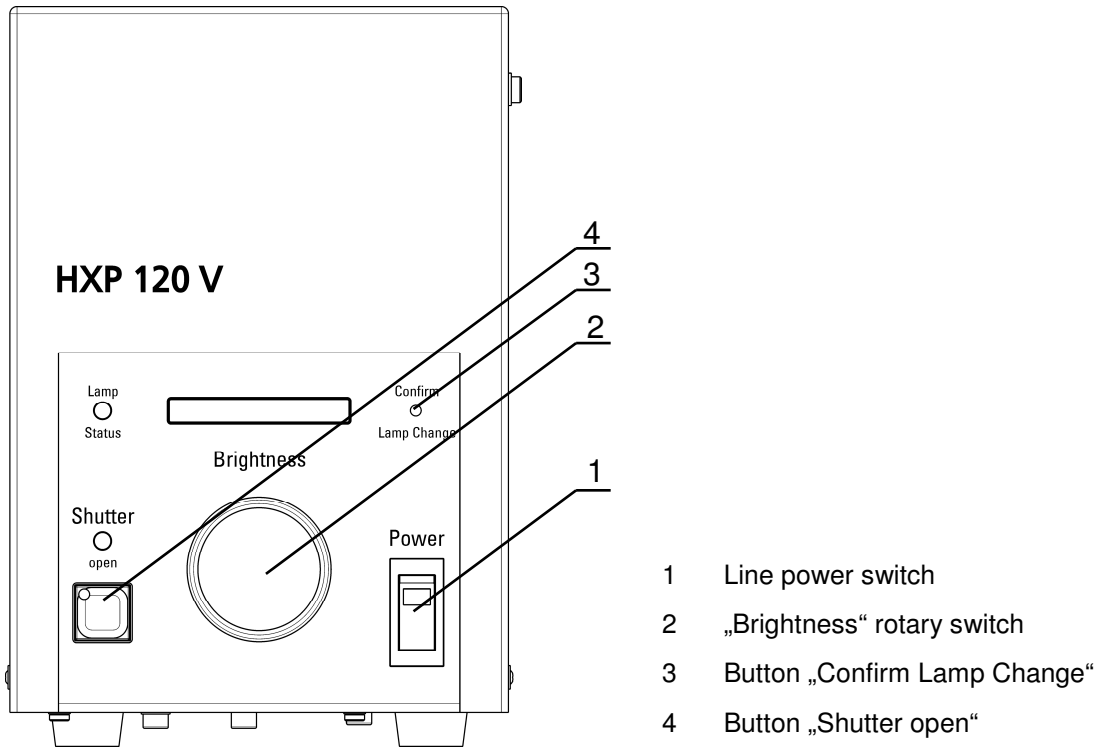
Figure 6: View of controls of the Compact Light Source HXP 120 V


The device is switched on and off at the main switch (Figure 6[1]). When the device is switched on the initialization runs. Both the dimmer unit and the integrated shutter move during this process.

The brightness emitted by the light source can be set to 20 different steps at the „Brightness“ rotary switch (Fig. 6[2]). In addition to the visual display of the brightness setting by the LED bar graph an audible signal is sounded at both extreme positions.

Pressing and holding (> 2sec) the „Confirm Lamp Change“ button (Fig. 6[3]) resets the internal lamp operating hours counter of the device. The reset is acknowledged by a beep. The lamp operating hours counter should be reset after every lamp change. The built-in lamp must in no case be operated longer than the maximum life specified by the manufacturer (see the lamp manufacturer’s instructions).

The „Shutter open“ button (Fig. 6[4]) controls the integrated shutter.

5.2 Indicators

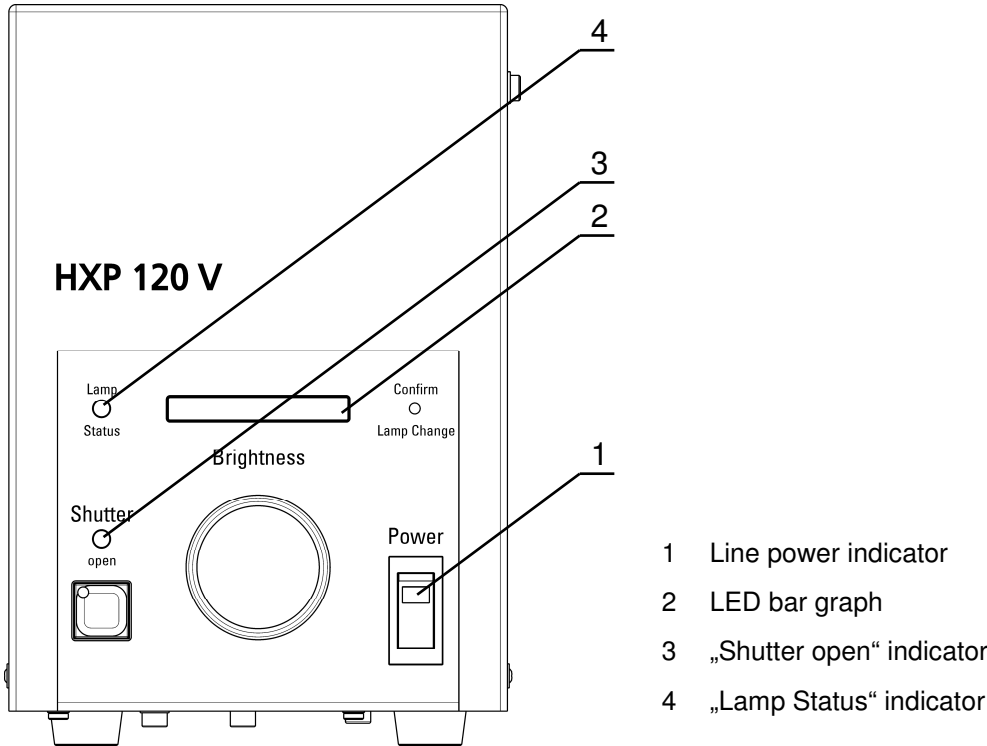- 1 Line power indicator
- 2 LED bar graph
- 3 „Shutter open“ indicator
- 4 „Lamp Status“ indicator

Figure 7: View of the indicators of the Compact Light Source HXP 120 V

The compact light source HXP 120 V has the following indicators:

- Power indicator (in the mains power switch, green, Fig. 7[1])
Lights when the device is switched on
- LED bar graph indicator (Fig. 7[2])
This graph indicates the brightness step emitted by the device. The longer the displayed LED bar, the higher the brightness is set.
- „Shutter open“ indicator (LED blue, Fig. 7[3])
Lights when the shutter integrated in the device is open.
- „Lamp Status“ indicator (LED multi-colour, Fig. 7[4])
After the lamp has fired, this LED indicates the state of the internal operation hours counter and therefore the age of the lamp in different colours.

„Lamp Status“ display	Internal operation hours counter	Comment
Green	< 2000 h	
Yellow	>2000 h	Prepare lamp change
Red	>2500 h	Change lamp
Flashing red		Lamp defective

Note: When the „Lamp Status“ display signals a defective lamp (flashing red), the device can be operated again after it has been shut off at the power switch (> 5sec) and the lamp operation hours counter is reset (by pressing the „Confirm Lamp Change“ button (Fig. 6[3])) after the device is switched on again.

- „Dimming end position“ indication (audible)
If the dimming is set to either end position by turning the „Brightness“ rotary switch (Fig. 6[2]) in, an audible signal (beep) is sounded.

5.3 Troubleshooting and remedies

Table 4 lists faults and defects which can occur during operation of the Compact Light Source HXP 120 V. The table also suggests procedures to remedy faults.

Problem	Possible defect	Remedy
Power indicator (Fig. 7[1]) does not light when the device is switched on, device shows no function	No power supply; power cable not connected correctly or defective	Inspect mains power and power cable
	Fuses in the device (Fig. 5[6]) defective	Replace fuses, use only approved fuse type
Power indicator (Fig. 7[1]) lights when the device is switched on, device shows no function	Device cover loose or removed	Firmly install / close device cover
	No lamp installed	Install lamp
Power indicator (Fig. 7[1]) lights when the device is switched on, „Lamp Status“ indicator (Fig. 7[4]) flashes red	Lamp life at end	Replace lamp and reset lamp operation hours counter (press "Confirm Lamp Change" button (Fig. 6[3]))
No / not enough light reaches the target system	Shutter closed ("Shutter open" indicator off) or „Brightness“ rotary switch at Min position	Open shutter at the appropriate switch (Fig. 6[2/4])
	Lamp not installed correctly	Check
	Optical fiber not properly plugged in the light output	Check
	Optical fiber defective / damaged	Replace / repair


Table 4: Faults and fault remedies

If the defect or fault cannot be remedied by applying the suggestions in Table 4, return the compact light source for service by the manufacturer giving a detailed description of the problem.

6 Service

6.1 Maintenance and cleaning

The Compact Light Source HXP 120 V does not need regular maintenance. This does not release the user from the obligation to carry out the statutory inspections for electric safety and accident prevention.

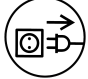 Switch off the Compact Light Source HXP 120 V before cleaning it. Use a dry, lintfree cloth for cleaning. When cleaning the device, avoid the entry of dust or other material through the ventilation slots. If the device is damaged (including external damage), switch it off immediately and return it to service.

If the Compact Light Source HXP 120 V is not used for a longer time, cover it suitably (e.g., use the plastic bag from the original packaging) to prevent the entry of dust.

6.2 Spare parts and accessories

 The user of the Compact Light Source HXP 120 V should only change the device fuses, the lamp and the glass of the protective filter. Removal and installation of the lamp and the protective filter are described in detail in chapters „4.2 Installing and changing the lamp“ and „4.3 Protective filter“. Sources from which to obtain the filter glass can be inquired from the supplier of the Compact Light Source HXP 120 V.

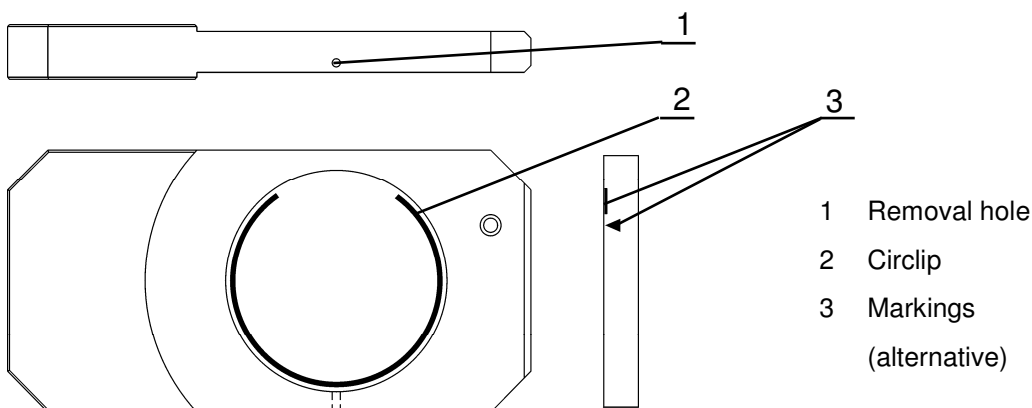
Figure 8: Protective filter and filter glass

Do not touch the filter glass with bare hands, sweat and fat will burn in the glass and render it unusable. The filter glass of the protective filter can be removed when you straighten a paper clip or use another suitable length of wire, insert it through the removal hole (Figure 8[1]) of the protective filter and compress the circlip (Figure 8[2]). The new filter glass can be installed in the filter after the old filter glass is removed. Make sure that the mark on the edge of the filter glass (Figure 8[3]) is oriented towards the lamp after the filter glass is installed. Installation of the circlip (Figure 8[2]) (when installed, the opening of the circlip should be opposite the removal hole) secures the filter glass in the protective filter and can be installed in the device. The loose fit of the filter glass is intentional.

To change the fuses, disconnect the connecting cable from the Compact Light Source HXP 120 V. The fuses are installed in a drawer at the device's line power input. Open it with a narrow screwdriver blade. Replace the fuses. Only the following fuse types are permitted as replacement:


Device fuse 5 x 20 2,5A, slow, switching capacity H	Manufacturer	Internet address
Series 181	Wickmann	www.wickmann.com
alternatively:		
Series SPT 5x20	Schurter	www.schurter.ch

Table 5: Permitted fuses for use with the Compact Light Source HXP 120 V

6.3 Repairs

No repair or maintenance except as specified in these operating instructions (lamp change; filter glass change; fuse change) must be carried out on the device. Repairs must only be carried out by the device supplier or an authorized repair center.

If the Compact Light Source HXP 120 V is defective and must be returned to the supplier or authorized repair center, use the original packaging of the device.

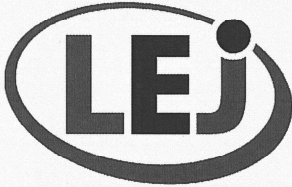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We declare the compliance of the device with the requirements of the Council Directive 2006 / 95 / EC - Low Voltage Directive and with the Council Directive 2004 / 108 / EC about the Electromagnetic Compatibility. Any modification of the product, not authorized by us, will invalidate this declaration.

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Leistungselektronik JENA GmbH